# PHOENIX ELECTRODE COMPANY PERCHLORATE ION ELECTRODES INSTRUCTION MANUAL

#### GENERAL INSTRUCTIONS

#### Introduction

The pHoenix Perchlorate Ion Electrodes are used to quickly, simply, accurately, and economically measure perchlorate in aqueous solutions.

#### Required Equipment

- 1
- A pH/mV meter or an ion meter, either line operated or portable.
- 2. Semi-logarithmic 4-cycle graph paper for preparing c alibration curves when using the meter in the mV mode.
- 3. A magnetic stirrer.
- 4.

The pHoenix Perchlorate Ion Electrode, Cat. No. PER1501, (reference electrode necessary), the pHoenix Perchlorate Ion Combination Glass Electrode, Cat. No. PER1502, or the Perchlorate Ion Combination Epoxy Electrode, Cat. No. PER1503.

The pHoenix Double Junction Reference Electrode, Cat. No. 5731429 (for use with the PER1501) with pHoenix Reference Filling Solution, Cat No. R001011, in the inner junction and with pHoenix Reference Filling Solution, Cat. No. R001044, in the outer junction.

#### Required Solutions

- 1. Deionized or distilled water for solution and standard p reparation.
- 2. pHoenix Ionic Strength Adjuster (ISA),  $2M (NH_4)_2SO_4$ , Cat. No. PERISO1. To prepare this solution from your own laborator y stock, half fill a 1000 ml volumetric flask with distilled w ater and add 264 grams of reagent-grade ammonium sulfate, ( $NH_4$ ) $_2SO_4$ . Swirl the flask gently to dissolve the solid. Fill the flask to the mark with distilled water, cap, and upend severa l times to mix the contents. ISA is added at the rate of 2 ml

of ISA to each 100 ml of standard or sample to adjust the ion ic strength to about 0.12M.

- 3. pHoenix Electrode Filling Solution,  $0.1M~(NH_4)_2SO_4/0.01M~KCl$ , Cat. No. R001045, for the PER1503 epoxy combination elect rode, or pHoenix Reference Filling Solution,  $0.1M~(NH_4)_2SO_4$ , Cat. No. R001044, for the 5731429 double junction electrode or PER1502 glass combination electrode.
- pHoenix Perchlorate Standard, 0.1M NaClO<sub>4</sub>, Cat. No. P ERASO1. To prepare this solution from your own laboratory sto ck, pipet 200 ml standard 0.5 M sodium hydroxide into a 600 ml beaker. Titrate to pH 7.00 against 60% perchloric acid using a calibrated pH electrode as the endpoint detector. Transfer, with rinsings, to a one liter volumetric flask and dilute to volume with distilled water. NOTE: Perchloric acid is explosive in contact with reducible substances and is a strong oxidizing agent. Use extreme caution when handling.
- 5. pHoenix Perchlorate Standard, 1000 ppm  ${\rm ClO_4}^{-1}$ , Cat. No.PER AS02. To prepare this solution from your own laboratory stock , pipet 20 ml standard 0.5 M sodium hydroxide into 200 ml dist illed water in a 600 ml beaker. Titrate to pH 7.00 against 60 % perchloric acid using a calibrated pH electrode as the endpo int indicator. Transfer, with rinsings, to a one liter volume tric flask and dilute to volume with distilled water. **NOTE**: P erchloric acid is explosive in contact with reducible substanc es and is a strong oxidizing agent. Use extreme caution when handling.

#### GENERAL PREPARATION

#### Electrode Preparation

Remove the rubber caps covering the electrode tips and any rubber i nserts covering the filling holes of the reference electrode. Fill the combination electrode or the reference electrode with the fill ing solution shipped with the electrode to a level just below the fill hole. No preparation is required for a sealed reference electrode. Gently shake the electrode downward in the same manner as a clinical thermometer to remove any air bubbles which may be trapped behind the perchlorate membrane. Prior to first usage, or after long-term storage, immerse the perchlorate membrane in perchlorate st andard for thirty minutes. The electrode is now ready for use.

Connect the electrodes to the proper terminals of the meter as recommended by the meter manufacturer.

#### 

- 1.
- To a 150 ml beaker, add 100 ml of distilled water and tw o ml of ISA. Place the beaker on a magnetic stirrer and begin stirring at a constant rate. After assuring that the meter is in the millivolt mode, lower the electrode tips into the solution.
- 2.

Using a pipet, add 1 ml of 0.1M or 1000 ppm perchlorate standard to the beaker. When the reading has stabilized, record the mV reading.

3.

Using a pipet, add 10 ml of the same perchlorate standard used above to the beaker. When the reading has stabilized, record the mV reading.

4

Determine the difference between the two readings. The e lectrode is operating correctly if the mV potential has change d by 56 " 2 mV assuming the solution temperature is between 20  $^{\circ}$  and 25 $^{\circ}$ C. See the **TROUBLESHOOTING** section if the potential c hange is not within this range.

**Slope** is defined as the change in potential observed when the concentration changes by a factor of 10.

# Electrode Slope Check (with ion meter) (check electrodes each day)

- 1.
- Prepare standard perchlorate solutions whose concentrations vary by ten fold. Use either the 0.1M or 1000 ppm perchlorate standard. Use the serial dilution method for this preparation.
- 2.

To a 150 ml beaker, add 100 ml of the lower value standard and 2 ml of ISA. Place the beaker on the magnetic stirrer and begin stirring at a constant rate. Lower the electrode tips into the solution. Assure that the meter is in the concentration mode.

3.

Adjust the meter to the concentration of the standard an d fix the value in the memory according to the meter manufactu rer's instructions.

- 4. Rinse the electrodes and blot dry.
- 5.

To another 150 ml beaker, add 100 ml of the higher value standard and 2 ml of ISA. Place the beaker on the magnetic s tirrer and begin stirring at a constant rate. Lower the elect rode tips into the solution.

6.

Adjust the meter to the concentration of the standard an d fix the value in the memory.

7.

Read the electrode slope according to the meter manufacturer's instructions. Correct electrode operation is in dicated by a slope of 90-100%. See the **TROUBLESHOOTING** sections if the slope is not within this range.

#### **MEASUREMENT**

#### Measuring Hints

All samples and standards should be at the same temperature for precise measurement. A difference of  $1^{\circ}\text{C}$  in temperature will result in a 2% measurement error.

The sensing membrane is normally subject to water uptake and might appear milky. This has no effect on performance.

Constant, but not violent, stirring is necessary for accurate measu rement. Magnetic stirrers can generate sufficient heat to change the solution temperature. To counteract this effect, place a piece of insulating material, such as styrofoam sheet, between the stirrer and beaker.

Always rinse the electrodes with distilled water and blot dry betwe en measurements. Use a clean, dry tissue to prevent cross-contamination.

For samples with high ionic strength, prepare standards whose composition is similar to the sample.

Always check to see that the membrane is free from air bubbles afte r immersion into standard or sample.

A slow responding electrode may be caused by interferences to the e lectrode. To restore proper performance, soak the electrode in dis tilled water for about 5 minutes to clean the membrane, rinse, and soak in standard solution for about 5 minutes.

#### Sample Requirements

All samples must be aqueous and not contain organics which can diss olve in the membrane or extract out the liquid ion exchanger.

The temperature of the standard solutions and of the sample solutions should be the same and below  $40^{\circ}$ C. About a 2% error will be introduced for a 1°C difference in temperature.

Interferences should be absent. If they are present, use the proce dures found in the **Interferences** section to remove them.

The pH range for the perchlorate ion electrode is 2.5-11. Neutrali ze samples outside this range with acid or base to bring them in range.

#### Units of Measurement

Perchlorate concentrations are measured in units of ppm as sodium p erchlorate, ppm as perchlorate, moles per liter, or any other conve nient concentration unit. Table 1 indicates some of the concentration units.

TABLE 1: Concentration Unit Conversion Factors

moles/liter	$\underline{\text{ppm ClO}_4}^{-1}$	$\underline{\text{ppm NaClO}_4}$
1.0X10 <sup>-2</sup>	995.0	1225.0
1.0X10 <sup>-3</sup>	99.5	122.5
1.0X10 <sup>-4</sup>	10.0	12.2

#### MEASUREMENT PROCEDURE

#### Direct Measurement

Direct measurement is a simple procedure for measuring a large numb er of samples. A single meter reading is all that is required for each sample. The ionic strength of samples and standards should be made the same by adjustment with ISA for all perchlorate solutions. The temperature of both sample solution and standard solutions s hould be the same.

#### Direct Measurement of Perchlorate (using a pH/mV meter)

1

By serial dilution, prepare three standard solutions from the 0.1 M or 1000 ppm stock standard. The resultant concentration should be  $10^{-2}$ ,  $10^{-3}$ , and  $10^{-4}\mathrm{M}$  or 1000, 100 and 10 ppm perchlorate standards. Add 2 ml of ISA per 100 ml of standard. Prepare standards with a composition similar to the samples if the samples have an ionic strength above 0.1M.

2.

Place the most dilute solution (10<sup>-4</sup>M or 10 ppm) in a 150 ml beaker on the magnetic stirrer and begin stirring at a con stant rate. After assuring that the meter is in the mV mode, lower the electrode tips into the solution. After the reading has stabilized, record the mV reading.

3.

Place the mid-range solution  $(10^{-3} \text{M or } 100 \text{ ppm})$  in a 150 ml beaker on the magnetic stirrer and begin stirring. After r insing the electrodes in distilled water, blot dry and immerse the electrode tips in the solution. When the reading has stabilized, record the mV value.

4.

Place the most concentrated solution  $(10^{-2}\text{M} \text{ or } 1000 \text{ ppm})$  in a 150 ml beaker on the magnetic stirrer and begin stirring. After rinsing the electrodes in distilled water, blot dry, a nd immerse the electrode tips in the solution. When the reading has stabilized, record the mV reading.

5.

Using the semi-logarithmic graph paper, plot the mV reading (linear axis) against the concentration (log axis). A typical calibration curve can be found in Figure 1. Extrapolate the calibration curve down to about  $1.0 \times 10^{-5} \mathrm{M}$  ( $1.0 \mathrm{~ppm~ClO_4^{-1}}$ ).

A calibration curve is constructed on semi-logarithmic paper when u sing the pH/mV meter in the millivolt mode. The measured elec trode potential in mV (linear axis) is plotted against the standard concentration (log axis). In the linear region of the curve, only three standards are necessary to determine a calibration curve. In the non-linear region, additional points must be measured. The direct measurement procedures given are for the linear portion of the curve. The non-linear portion of the curve requires the use of low level procedures.

б.

To a clean, dry 150 ml beaker, add 100 ml of the sample and 2 ml of ISA. Place the beaker on the magnetic stirrer and begin stirring at a constant rate. Rinse the electrode tips with distilled water, blot dry, and lower the electrode tips into the solution. When the reading has stabilized, record the mV reading. Using the calibration curve, determine the sample concentration.

7.

The calibration should be checked every two hours. Assu ming no change in ambient temperature, place the electrode tip s in the mid-range standard. After the reading has stabilized, compare it to the original reading recorded in Step 3 above. A reading differing by more than 0.5 mV or a change in the a mbient temperature will necessitate the repetition of Steps 2-5 above. A new calibration curve should be prepared daily.

#### Direct Measurement of Perchlorate (using an ion meter)

1.

By serial dilution of the 0.1M or 1000 ppm perchlorate s tandard, prepare two perchlorate standards whose concentration is near the expected sample concentration. Measure out 100 m l of each standard into individual 150 ml beakers and add 2 ml of ISA to each.

2. .

Place the most dilute solution on the magnetic stirrer a nd begin stirring at a constant rate. Assure that the meter i s in the concentration mode. Lower the electrode tips into the solution.

3

Adjust the meter to the concentration of the perchlorate standard and fix the value in the memory according to the meter manufacturer's instructions after stabilization of the reading.

4. Rinse the electrodes with distilled water and blot dry.

- Place the more concentrated solution on the magnetic stirrer and begin stirring at a constant rate. Lower the electrode tips into the solution.
- Adjust the meter to the concentration of the perchlorate standard and fix the value in the memory according to the meter manufacturer's instructions after stabilization of the reading.
- For low level measurements, place the rinsed, dried elec trodes into a solution containing 100 ml of distilled water and 2 ml of ISA. After stabilization, fix the blank value in the meter according to the meter manufacturer's instructions.
- Place 100 ml of the sample and 2 ml of ISA in a 150 ml b eaker. Place the beaker on the magnetic stirrer and begin stirring.
- Immerse the electrode tips in the solution and wait for the reading to stabilize. Read the concentration directly from the meter display.
- The calibration should be checked every two hours. Assu ming no change in ambient temperature, place the electrode tip s in the first perchlorate standard. After the reading has s tabilized, compare it to the original reading in Step 3 above. A reading differing by more than 0.5 mV or a change in ambie nt temperature will necessitate the repetition of Steps 2-6 ab ove. The meter should be re-calibrated daily.

#### Low Level Perchlorate Determination (using a pH/mV meter)

This procedure is recommended for solutions with ionic strengths le ss than  $1.0 \times 10^{-2} \text{M}$ . If the solution is high in ionic strength, but l ow in perchlorate, use the same procedure, but prepare a calibration solution with a composition similar to the sample.

- 1. Using 20 ml of standard ISA, dilute to 100 ml with distilled water. This low level ISA  $[0.4M\ (NH_4)_2SO_4]$  is added at the rate of 1 ml low level ISA to each 100 ml of sample. The background ionic strength will be  $4.0\times10^{-3}M$ .
- 2. Dilute 1 ml of 0.1M standard to 100 ml to prepare a  $1.0 \rm X 10^{-3} M \ ClO_4^{-1}$  solution for measurements in moles per liter.

Use the 1000 ppm standard for preparing a 100 ppm  ${\rm ClO_4}^{-1}$  standard by diluting 10 ml of the 1000 ppm standard to 100 ml. Standards should be prepared fresh daily.

3.

Add 1 ml of the low level ISA to a 100 ml volumetric f lask and fill to the mark with distilled water. Pour this solution into a 150 ml beaker and place the beaker on the magnetic stirrer. Begin stirring at a constant rate.

4.

Place the electrode tips in the solution. Assure that the meter is in the mV mode.

5.

Add increments of the  $1.0 \times 10^{-3} \rm M$  or 100 ppm standards as g iven in Table 2 below.

6.

After the reading has stabilized, record the mV reading after each addition.

TABLE 2: Step-wise Calibration for Low Level Perchlorate
Measurements

	Added			tration
Step	<u>Pipet</u>	Volume (ml)	$\underline{M}\ \mathtt{Clo}_{\underline{4}}^{-1}$	$\underline{\text{ppm ClO}_4}^{-1}$
1	A	0.1	1.0X10 <sup>-6</sup>	0.1
2	A	0.1	$2.0 \times 10^{-6}$	0.2
3	A	0.2	$4.0 \times 10^{-6}$	0.4
4	A	0.2	$6.0 \times 10^{-6}$	0.6
5	A	0.4	$9.9 \times 10^{-6}$	1.0
6	В	2.0	$2.9 \times 10^{-5}$	2.9
7	В	2.0	4.8X10 <sup>-5</sup>	4.8

Pipet A = 1 ml graduated pipet

Pipet B = 2 ml pipet

Solutions: additions of 1.0X10<sup>-3</sup>M or 100 ppm standard to 100 ml of ISA as prepared in Step 3 above.

7.

On semi-logarithmic graph paper, plot the millivolt  $\,$  r eading (linear axis) against the concentration (log axis) as i  $\,$  n Figure 1.

- 8. Rinse the electrodes with distilled water and blot dry.
- 9. Measure out 100 ml of the sample into a 150 ml beaker, a

dd 1 ml of low level ISA. Place the beaker on the magnetic st irrer and begin stirring. Lower the electrode tips into the s olution. After the reading has stabilized, record the mV reading and determine the concentration from the low level calibration curve.

10.

Prepare a new low level calibration curve daily. Check the calibration curve every two hours by repeating Steps 2-7.

#### Low Level Perchlorate Determination (using an ion meter)

Follow the procedure given for normal perchlorate determinations us ing an ion meter and the blank correction procedure.

#### ELECTRODE CHARACTERISTICS

#### Reproducibility

Electrode measurements reproducible to "2% can be obtained if the e lectrode is calibrated every hour. Factors such as temperature flu ctuations, drift, and noise limit reproducibility. Reproducibility is independent of concentration within the electrode's operating range.

#### Interferences

Certain anions are electrode interferences and will cause electrode malfunction, drift or measurement errors if present in high enough levels. The level of interfering common anions that will cause a 10% error at three levels of perchlorate is given in Table 3.

TABLE 3: Concentration of Possible Interferences Causing a 10% Error at Various Levels of Perchlorate; Background Ionic Strength of 0.12M (NH<sub>4</sub>)<sub>2</sub>SO<sub>4</sub>.

Interference (moles/liter		$\underline{10^{-3}M}$	$10^{-4} \text{M ClO}_4^{-1}$
$NO_3^{-1}$	$5.0 \times 10^{-1}$	$5.0 \times 10^{-2}$	$5.0 \times 10^{-3}$
$\mathrm{NO_2}^{-1}$	$5.0 \times 10^{-1}$	$1.0 \times 10^{-2}$	$1.0 \times 10^{-3}$
$\mathtt{Br}^{-1}$	$4.0 \times 10^{-1}$	$4.0 \times 10^{-2}$	$4.0 \times 10^{-3}$
$CN^{-1}$	$4.0 \times 10^{-1}$	$4.0 \times 10^{-2}$	$4.0 \times 10^{-3}$
${\rm ClO_3}^{-1}$	$2.0x10^{-1}$	$2.0 \times 10^{-2}$	$2.0 \times 10^{-3}$
$\mathtt{I}^{-1}$	$2.0 \times 10^{-2}$	$2.0 \times 10^{-3}$	$2.0 \times 10^{-4}$

Interferences	5		
(ppm) NO <sub>3</sub> -1	$\frac{100 \text{ ppm}}{3.1 \times 10^3}$	$\frac{10 \text{ ppm}}{3.1 \text{x} 10^2}$	$\frac{1 \text{ ppm}}{3.1 \text{x} 10^{1}}$
$NO_2^{-1}$	$2.3x10^{3}$	$2.3x10^{2}$	$2.3x10^{1}$
$\mathrm{Br}^{-1}$	$3.2x10^{3}$	$3.2x10^{2}$	$3.2x10^{1}$
$CN^{-1}$	$1.0 \times 10^{3}$	$1.0 \times 10^{2}$	$1.2 \times 10^{1}$
${\rm ClO_3}^{-1}$	$1.7 \times 10^3$	$1.7x10^{2}$	$1.7x10^{1}$
$\mathtt{I}^{-1}$	$2.5x10^{2}$	$2.5 \text{x} 10^{1}$	2.5

Interferences such as bromide, iodide and cyanide can be removed by precipitation with 0.5 grams of silver sulfate added to 100 ml of sample. Nitrite interference can be removed by adding 0.3 grams of sulfamic acid to 100 ml of sample.

The above interference removal procedures require similar treatment of standards as well as samples.

If the electrode is exposed to high levels of interfering ions which cannot be removed, the electrode reading may drift and the

response may become sluggish. Restore performance by soaking in distilled water for 30 minutes followed by soaking in perchlorate standard for 30 minutes.

#### Temperature Influences

Samples and standards should be at the same temperature, since elec trode potentials are influenced by changes in temperature. A  $1^{\circ}\text{C}$  difference in temperature results in a 2% error at the  $10^{-3}\text{M}$  level. B ecause of the solubility equilibria on which the electrode depends, the absolute potential of the reference electrode changes slowly w ith temperature. The slope of the perchlorate electrode, as indicated by the factor "S" in the Nernst equation, also varies with temperature. Table 4 gives values for the "S" factor in the Nernst equation for the perchlorate ion.

The operating range of the perchlorate ion electrode is  $0^{\circ}-40^{\circ}\text{C}$ , provided that temperature equilibrium has occurred.

If the temperature varies substantially from room temperature, equilibrium times up to one hour are recommended.

TABLE 4: Temperature vs. Values for the Electrode Slope

Temp ( $^{\circ}$ C) "S"

0	54.20
10	56.18
20	58.18
25	59.16
30	60.15
40	62.13
50	64.11

#### Electrode Response

Plotting the mV potential against the perchlorate concentration on semi-logarithmic paper results in a straight line with a slope of a bout 56 mV per decade. (Refer to Figure 1.)

The time needed to reach 99% of the stable electrode potential read ing, the electrode response time, varies from one minute or less in highly concentrated solutions to several minutes near the detection limit. (Refer to Figure 2.)

#### Limits of Detection

The upper limit of detection in pure sodium perchlorate solutions is 1M. In the presence of other ions, the upper limit of detection is above  $10^{-1}M$  perchlorate, but two factors influence this upper limit. Both the possibility of a liquid junction potential developing at the reference electrode and the salt extraction effect influence this upper limit. Some salts may infuse into the electrode membrane at high salt concentrations, causing deviation from the theoretical response. Either dilute samples between 1M and  $10^{-1}M$  or calibrate the electrode at 4 or 5 intermediate points.

The lower limit of detection is influenced by the slight water solu bility of the ion exchanger used in the sensing portion of the elec trode. Refer to Figure 1 for a comparison of the theoretical response to the actual response at low levels of perchlorate. Perchlora te measurements below  $10^{-5} \rm M~ClO_4^{-1}~(1.0~ppm~as~ClO_4^{-1})$  should employ low level procedures.

#### pH Effects

The operating range of the perchlorate electrode is from pH 2.5 to pH 11.

#### Electrode Life

The perchlorate electrode will last six months in normal laboratory use. On-line measurement might shorten operational lifetime to se veral months. In time, the response time will increase and the cal ibration slope will decrease to the point calibration is difficult and electrode replacement is required.

#### Electrode Storage

The perchlorate electrodes may be stored for short periods of time in  $10^{-2}$ M perchlorate solution. For longer storage (longer than two weeks), rinse and dry the perchlorate membrane and cover the tip with any protective cap shipped with the electrodes. The reference portion of the combination electrode (or the outer chamber of the reference electrode) should be drained of filling solution, if refill able and the rubber insert placed over the filling hole.

#### ELECTRODE THEORY

#### Electrode Operation

The perchlorate electrode consists of an electrode body containing a liquid internal filling solution in contact with a gelled organop hilic membrane containing a perchlorate ion exchanger. When the membrane is in contact with a solution containing free perchlorate ions, an electrode potential develops across the membrane. This electrode potential is measured against a constant reference potential, using a standard pH/mV meter or an ion meter. The level of perchlorate ions, corresponding to the measured potential, is described by the Nernst equation:

 $E = E\$ - S \log X$ 

where: E = measured electrode potential

E\$= reference potential (a constant)

S = electrode slope (-56 mV/decade)

X = level of perchlorate ions in

solution

The activity, X, represents the effective concentration of the ions in solution. The total perchlorate ion concentration,  $C_t$ , is the s um of free perchlorate ion,  $C_f$ , and complexed or bound perchlorate ion,  $C_b$ . The electrode is able to respond to only the free ions, wh ose concentration is:

$$C_f = C_t - C_b$$

Since perchlorate ions form very few stable complexes, the free ion concentration may be equated to the total ion concentration.

The activity is related to the free ion concentration,  $C_f$ , by the a ctivity coefficient, a, by:

 $X = a C_f$ 

Activity coefficients vary, depending on total ions strength, I, de fined as:

 $I = 1/2 3C_xZ_x5$ 

where:  $C_x$  = concentration of ion X  $Z_x$  = charge of ion X 3 = sum of all of the types of ions in the solution

In the case of high and constant ionic strength relative to the sen sed ion concentration, the activity coefficient, a, is constant and the activity, X, is directly proportional to the concentration.

To adjust the background ionic strength to a high and constant value, ionic strength adjuster (ISA) is added to samples and standards. The recommended ISA for perchlorate is  $(NH_4)_2SO_4$ . Solutions other than this may be used as ionic strength adjusters as long as ions that they contain do not interfere with the electrode's response to perchlorate ions.

The reference electrode must also be considered. When two solution s of different composition are brought into contact with one anothe r, liquid junction potentials arise. Millivolt potentials occur fr om the inter-diffusion of ions in the two solutions. Electrode charge will be carried unequally across the solution boundary resulting in a potential difference between the two solutions, since ions diffuse at different rates. When making measurements, it is important to remember that this potential be the same when the reference is in the standardizing solution as well as in the sample solution or the change in liquid junction potential will appear as an error in the measured electrode potential.

The composition of the liquid junction filling solution in the reference electrode is most important. The speed with which the positive and negative ions in the filling solutions diffuse into the samples should be as nearly equal as possible, that is, the filling solution should be equitransferent. No junction potential can result if the rate at which positive and negative charge carried into the sample is equal.

Strongly acidic (pH = 0-2) and strongly basic (pH = 12-14) solution s are particularly troublesome to measure. The high mobility of hy drogen and hydroxide ions in samples make it impossible to mask the ir effect on the junction potential with any concentration of equit ransferent salt. One must either calibrate the electrodes in the s ame pH range as the sample or use a known incremental method for io n measurement.

#### TROUBLESHOOTING GUIDE

The goal of troubleshooting is the isolation of a problem through c hecking each of the system components in turn: the meter, the plast ic-ware, the electrodes, the standards & reagents, the sample, and the technique.

#### Meter

The meter may be checked by following the check-out procedure in the instrument instruction manual.

#### Plastic-ware

Clean plastic-ware is essential for good measurement. Be sure to w ash the plastic-ware well with a mild detergent and rinse very well with distilled or deionized water.

#### Electrodes

The electrodes may be checked by using the procedure found in the s ections entitled **Electrode Slope Check**.

- 1.

  Be sure to use distilled or deionized water when following the procedures given in **Electrode Slope Check**.
- 2.

  If the electrode fails to respond as expected, see the s ection **Measuring Hints**. Repeat the slope check.
- If the electrodes still fail to respond as expected, substitute another perchlorate ion electrode that is known to be in good working order for the questionable electrode. If the problem persists and you are using an electrode pair, t ry the same routine with a working reference electrode.
- 4.

  If the problem persists, the standards or reagent may be of poor quality, interferences in the sample may be present or the technique may be faulty. (See Standards & Reagents, S ample, and Technique sections below.)

5.

If another electrode is not available for test purposes, or if the electrode in use is suspect, review the instruction manual and be sure to:

- Clean and rinse the electrodes thoroughly.
- Prepare the electrodes properly.
- Use the proper filling solution.
- Adjust the pH and the ionic strength of the solution by use of the proper ISA.
- Measure correctly and accurately.
- Review TROUBLESHOOTING HINTS.

#### Standards & Reagents

Whenever problems arise with the measuring procedure that has been used successfully in the past, be sure to check the standard and re agent solutions. If in doubt about the credibility of any of the s olutions, prepare them again. Errors may result from contamination of the ISA, incorrect dilution of standards, poor quality distille d/deionized water, or a simple mathematical miscalculation.

#### Sample

Look for possible interferences, complexing agents, or substances we hich could affect the response or physically damage the sensing electrode (or the reference electrode) if the electrodes work perfectly in the standard, but not in the sample.

Try to determine the composition of the samples prior to testing to eliminate a problem before it starts. (See Measuring Hints, Sample Requirements, and Interferences.)

#### Technique

Be sure that the electrode's limit of detection has not been exceed ed. Be sure that the analysis method is clearly understood and is compatible with the sample. Refer to the instruction manual again. Reread **GENERAL PREPARATION** and **ELECTRODE CHARACTERISTICS**.

If trouble still persists, call pHoenix Electrode Company at 1-800-522-7920 and ask for the Technical Services Department.

### TROUBLESHOOTING HINTS

Symptom	Possible Causes	Next Step
Out of Range Reading		check meter with shorting strap see meter instruction
	defective electrode	ahadr alagtrada
		peration
plugged	electrodes not d in properly and	unplug electrodes reseat
		e be sure reference electrode is filled
	air bubble on membrane	remove bubble by re-dipping electrode
	electrodes not in solution	put electrodes in solution
Readings (reading continuously or	defective meter	check meter with shorting strap
	air bubble on membrane	remove bubble by re-dipping electrode
	defective electrode	replace electrode
	ISA not used	use recommended ISA
	meter or stirrer not grounded	ground meter or stirrer
	electrode exposed to interferences standard	soak electrode in perchlorate
	outer filling solution level too low	fill electrode to level just below the fill hole
Drift (reading slowly changing in one direction)	at different	dards allow solutions to come to room temper- ature before measurement

electrode exposed soak electrode in to interferences perchlorate

standard

incorrect reference use recommended filling solution filling solution

Low Slope or standards contaminprepare fresh No Sl

ated or incorrectly standards ope

made

ISA not used use recommended

ISA

standard used as ISA use ISA

electrode exposed soak electrode in to interferences perchlorate

standard

defective electrode check electrode operation

air bubble on remove bubble by membrane re-dipping probe

"Incorrect Answer" incorrect scaling plot millivolts on (but calibration of semi-log paper the linear axis. On cur ve is good) the log axis, be

sure concentration

numbers within each

decade are increa sing with increasing

concentration

incorrect sign be sure to note sign of millivolt number

correctly

incorrect standards prepare fresh standards

wrong units used apply correct conversion factor:

 $10^{-3}M = 122.5 \text{ ppm as}$ 

 $NaClO_4 = 99.5 ppm as$ 

 $ClO_{4}^{-1}$ 

sample carryover rinse electrodes thoroughly between

samples

#### **SPECIFICATIONS**

Concentration Range:  $1M \text{ to } 7 \times 10^{-6}M$ 

 $(1.0 \times 10^5 \text{ to } 0.7 \text{ ppm as } \text{ClO}_4^{-1})$ 

pH Range: 2.5 to 11

Temperature Range:  $0^{\circ}$  to  $40^{\circ}$ C

Resistance: 100 Mohms

Reproducibility: "2%

Samples: aqueous solutions only

no organic solvents

Size: 110 mm in length

12 mm in diameter 1 m cable length

Storage: Store in dilute perchlorate solution

#### ORDERING INFORMATION

P/N DESCRIPTION

PER1501 Perchlorate Electrode, mono (reference

electrode necessary), PVC body

PER1502 Perchlorate Electrode, combination, glass body

PER1503 Perchlorate Electrode, combination, epoxy body

5731429 Reference Electrode, double junction, sleeve, epoxy

body, for use with the PER1501

PERAS01 Perchlorate Standard, 0.1M NaClO<sub>4</sub>

PERAS02 Perchlorate Standard. 1000 ppm ClO<sub>4</sub><sup>-1</sup>

PERIS01 Perchlorate ISA (Ionic Strength Adjustor),

2M (NH<sub>4</sub>)<sub>2</sub>SO<sub>4</sub>

R001011 5731429 Reference Electrode Inner Filling Solution

4M KC1 (with Aq+)

R001044 5731429 Reference Electrode Outer Filling

Solution & PER1502 Glass Combination Electrode

Filling Solution,  $0.1M (NH_4)_2SO_4$ 

R001045 PER1503 Epoxy Combination Electrode Filling

Solution,  $0.1M (NH_4)_2SO_4/0.01M KCl (with Ag+)$ 

## TABLE OF CONTENTS

Gener	al Instructions 1 introduction 2 required equipment 2 required solutions 2	1 1
Gener	al Preparation	2
Measu	rement4measuring hints4sample requirements4units of measurement5	4 4
	direct measurement	5
	(using an ion meter)	9
Elect	rode Characteristics9reproducibility9interferences10temperature influences11electrode response12limits of detection12pH effects13electrode life13electrode storage13	9 0 1 2 2
Elect	rode Theory       electrode operation	
Trouk Speci Order	leshooting Guide	5 9